

Changing MCA bins in EXAFS code:

1. Exit EXAFS code using the STOP button (not the tiny stop-sign icon on the top).
2. Use the MCA utility to decide what the bins should be.
3. Edit the scaler-map file "2\_6602+2\_XIA.cfg" so that Scaler 2 has the correct values for roilow and roihigh.
4. Restart the EXAFS code by clicking on the white arrow (the one which goes black when it's running).
5. Reload the appropriate scan-parameter file.
6. Use the scan editor to make sure the regions are correct and the file names are what you want.